



PIXIS: 1300

The PIXIS: 1300 series of cameras from Princeton Instruments (PI) are fully integrated, low noise cameras designed expressly for quantitative scientific imaging applications. Choose front-illuminated (F), back-illuminated (B), eXcelon™ or back-illuminated deep depletion (BR) versions of the 1340 x 1300 CCD for optimized performance from UV to NIR. Designed utilizing PI's exclusive XP cooling technology, PIXIS are the only cameras that offer cooling up to -70°C, while the all-metal, hermetically sealed design, with the industry's only lifetime vacuum guarantee, ensures maintenance free operation. High QE and ultra low-noise electronics make the PIXIS: 1300 series of cameras ideal for demanding, low light applications such as astronomy, Bose-Einstein Condensate (BEC), solar cell inspection, chemiluminescence and fluorescence imaging. The eXcelon and back-illuminated, deep depletion technology of the PIXIS:1300 delivers the highest sensitivity in the NIR while suppressing the etaloning that occurs in standard back illuminated CCDs. Dual speed operation at 100kHz or 2MHz allow these cameras to be used for steady state as well as fast kinetics studies.

FEATURE	BENEFITS
eXcelon™ technology	Higher QE in the UV and near IR regions; Extremely low etaloning; 100x lower dark charge than back-illuminated deep depletion
1340 x 1300 imaging array; 20µm x 20µm pixels Scientific grade CCD	High spatial resolution and large field of view Low noise, few defects, linear response
All-metal, hermetic vacuum seals; Lifetime vacuum guarantee	No out-gassing (as in epoxy seals) which can compromise vacuum performance; Worry free operation
Deep cooling	Low dark noise allows detection of faint signals; CoolCUBE II, a compact room temperature coolant circulator, is available for vibration sensitive environments
Single fused silica vacuum window	Minimizes reflection losses from UV to IR; No optical losses due to multiple optical surfaces; Optional AR coating and wedge windows available
Optional UV phosphor coatings	Enhances sensitivity throughout the UV to below 200 nm (not available for eXcelon cameras)
Large 45 mm integrated shutter	No vignetting even at low f-numbers. Easily acquire bias and dark reference.
Low noise electronics	Best performance for low light level applications
Dual digitizers	Dual-speed digitization allows complete freedom to select between "slow operation" for low noise and highest SNR or "fast operation" for rapid image acquisition
Software selectable system gains	Flexibility to optimize signal-to-noise ratio and dynamic range
Kinetics	Custom readout mode offers microsecond resolution
USB2.0 data interface	Plug-and-play operation with desktops or laptops; Optional fiber optic interface for remote operation
LightField™ for 64-bit Windows 7 or WinSpec for 32-bit Windows	Flexible software packages for data acquisition, display and analysis; LightField offers intuitive, cutting edge user interface, IntelliCal™ and more
PICAM/PVCAM software development kits (SDKs)	Universal programming interfaces for easy custom programming; Compatible with Windows XP/Vista/7 and Linux

Applications:

semiconductor failure analysis, astronomy, photometry, laser beam profiling, luminescence and fluorescence imaging, and Bose-Einstein Condensate (BEC)

SPECIFICATIONS




	PIXIS: 1300F	PIXIS: 1300B_eXcelon	PIXIS: 1300B	PIXIS: 1300BR
Features	Front-illuminated CCD. Affordable technology for moderate light level applications. No etaloning.	Back-illuminated CCD. Highest sensitivity in the visible region. High sensitivity in the NIR. Extremely low etaloning. 100x lower dark charge than the BR.	Back-illuminated CCD. Highest sensitivity in the visible region.	Back-illuminated, deep depletion CCD. Ideal for NIR applications. Highest sensitivity and no etaloning.
CCD Image Sensor	Princeton Instruments' proprietary CCD, front-illuminated, grade 1, AIMO	Princeton Instruments' proprietary CCD, grade 1, AIMO	Princeton Instruments' proprietary CCD, back-illuminated, grade 1, AIMO	Princeton Instruments' proprietary CCD, back-illuminated deep depletion, grade 1, NIMO
Dark current @ -60°C (e-/p/sec)	0.01 (typical) 0.05 (max)	0.01 (typical) 0.05 (max)	0.01 (typical) 0.05 (max)	0.32 (typical) 0.65 (max)
CCD UV coating	Optional UV coating			
CCD format	1340 x 1300 imaging pixels; 20µm x 20µm pixels; 100% fill factor			
Imaging area	26.8 x 26 mm (optically centered)			
Lens mount	F-mount with integral 45mm shutter			
Deepest cooling temperature	< -70°C (typical), -60°C (guaranteed) with CoolCUBEII liquid circulator < -65°C (typical), -55°C (guaranteed) with air			
Thermostating precision	±0.05 °C			
Cooling method	Thermoelectric air or liquid cooling (CoolCUBE II required)			
Full well: Single pixel	250 ke- (typical), 200 ke- (min)			
Output node	1000 ke- (typical), 800 ke- (min)			
ADC speed/bits	100kHz/16-bit and 2MHz/16-bit			
System read noise @ 100kHz	2 e- rms (typical), 3 e- rms (max)			
@ 2MHz	12 e- rms (typical), 16 e- rms (max)			
Vertical shift speed	27 µsec/row (programmable)			
Non-linearity	<1% @ 100kHz			
Software selectable gains	1, 2, 4 e-/ADU (low noise output); 3.5, 7, 14 e-/ADU (high capacity output)			
Operating systems supported	Windows XP/Vista/7; Linux			
Data interface	USB2.0 (5m interface cable provided); Optional Fiberoptic interface is available for remote operation			
I/O signals	Two MCX connectors for programmable frame readout, shutter, trigger in			
Operating environment	+5 to +30°C non-condensing			
Certification	CE			
Dimensions / Weight	19.51cm (7.67") x 11.81cm (4.65") x 11.38cm (4.48") (L x W x H) / 2.5kg (5.5lbs)			

All specifications subject to change

FRAME RATE

	Readout Time	
	@ 2 MHz	@ 100 kHz
Binning 1 x 1	0.93 sec	16 sec
2 x 2	0.402 sec	4.762 sec
8 x 8	0.12 sec	0.491 sec
16 x 16	0.083 sec	0.209 sec


QE DATA

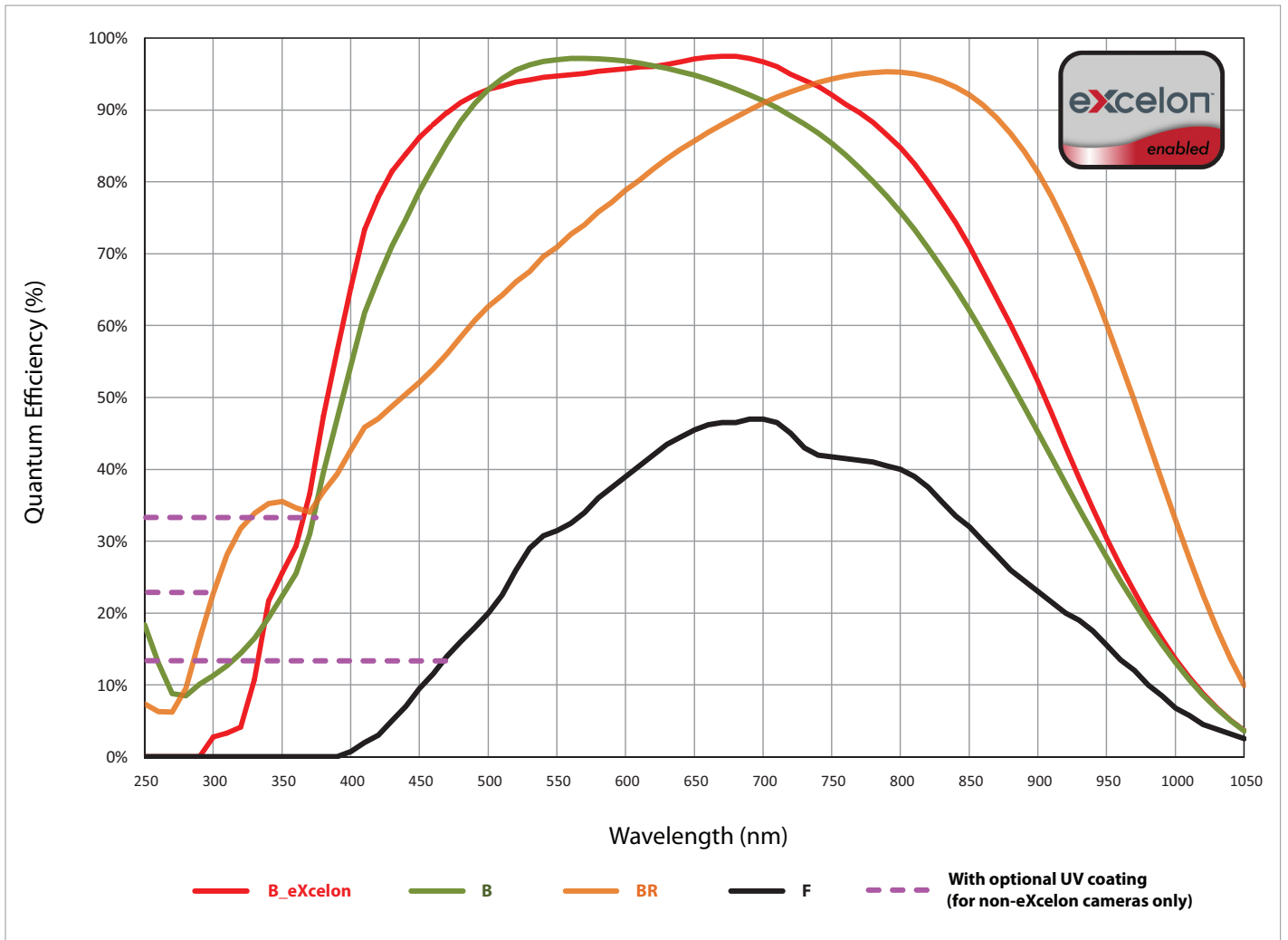

NOTE:

Graph shows typical Quantum Efficiency (QE) data measured at + 25°C. QE decreases at normal operating temperatures. For the best results for your application, please discuss the specific parameters of your experiment with your sales representative.


VACUUM WINDOW AR COATINGS

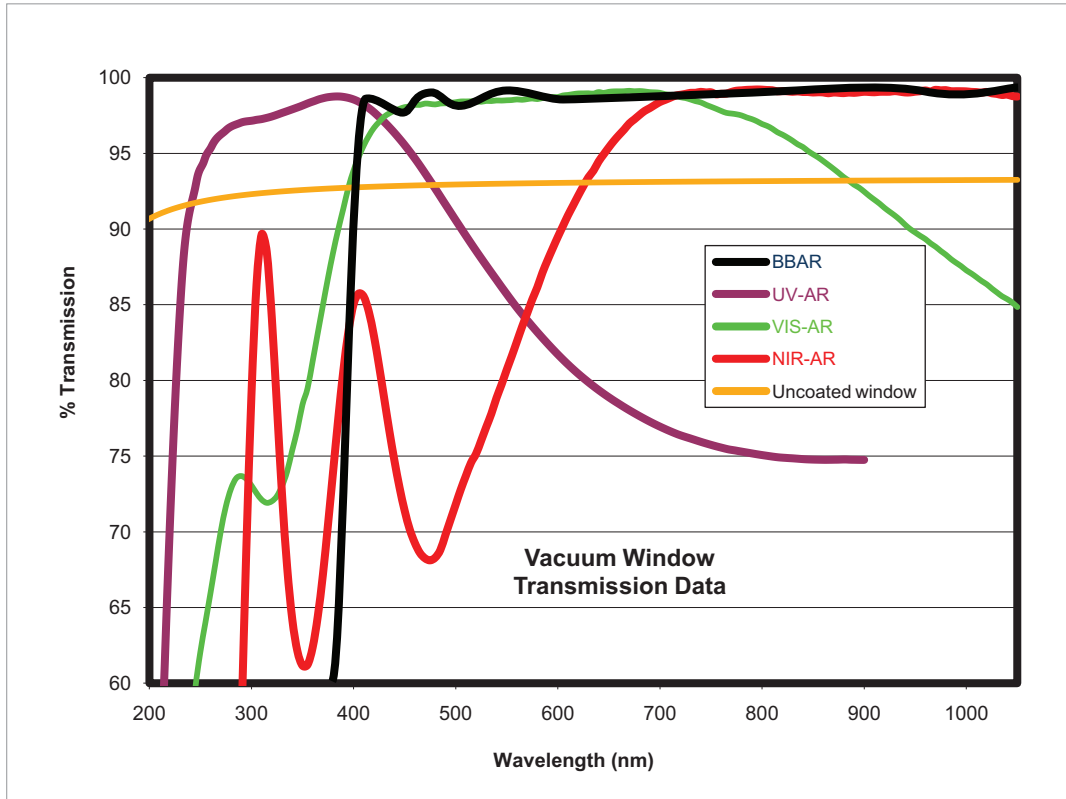


NOTE:

Standard anti-reflection (AR) coatings shown. Custom AR coatings and wedge window options are also available. Contact your local sales representative for more information.


eXcelon Performance

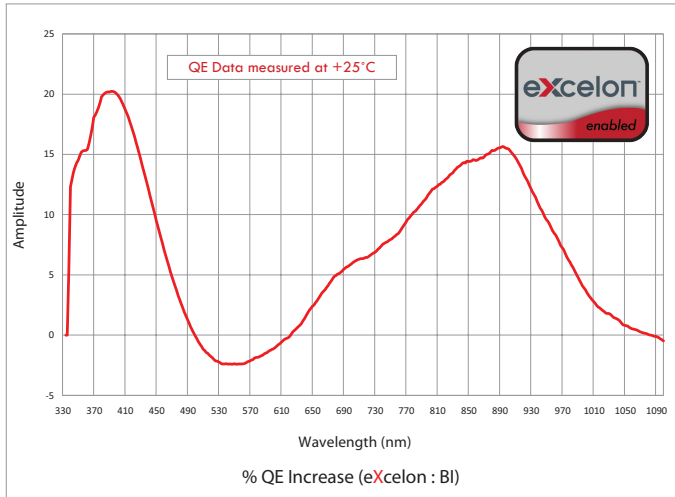

B_eXcelon provides superior QE over the standard back illuminated ("B") version in the UV-NIR range.

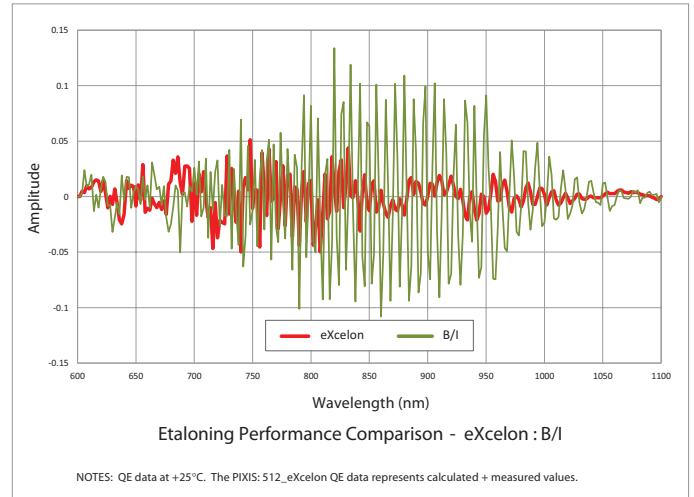


B_eXcelon provides significantly lower etaloning (unwanted fringes) compared to standard back illuminated ("B") version.

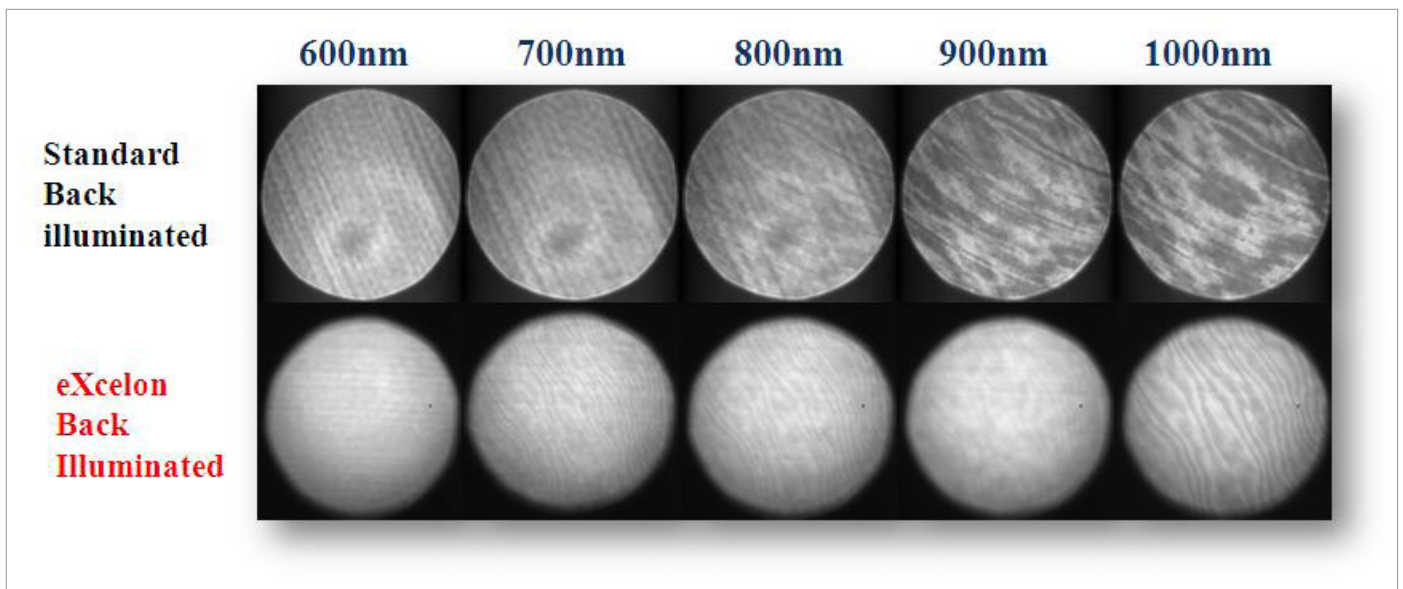


Data taken with white light source through a monochromator comparing etaloning performance of eXcelon vs. back-illuminated CCDs.

PIXIS: 1300 (AIR COOLED)

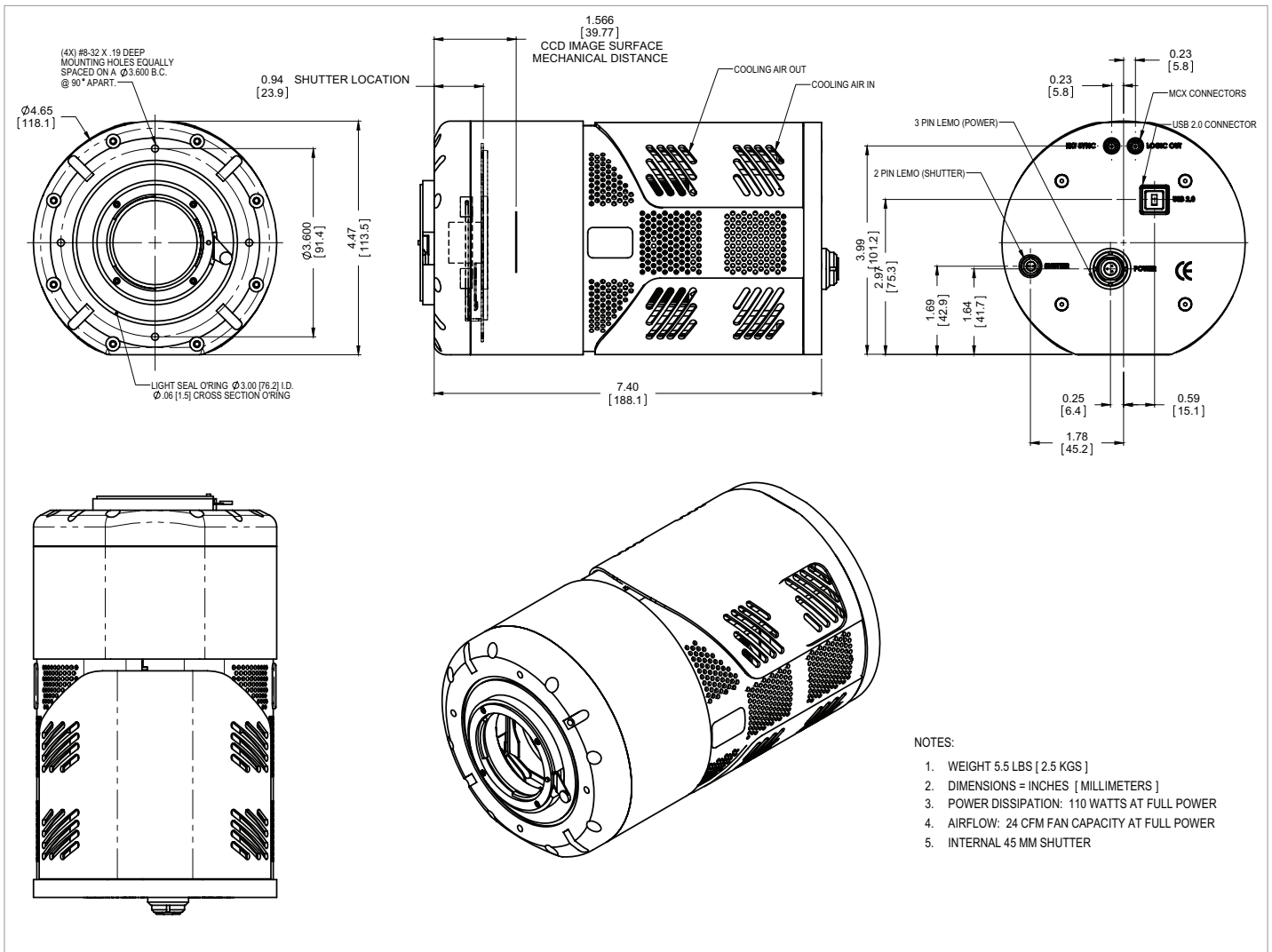

PIXIS: 1300 (LIQUID COOLED)

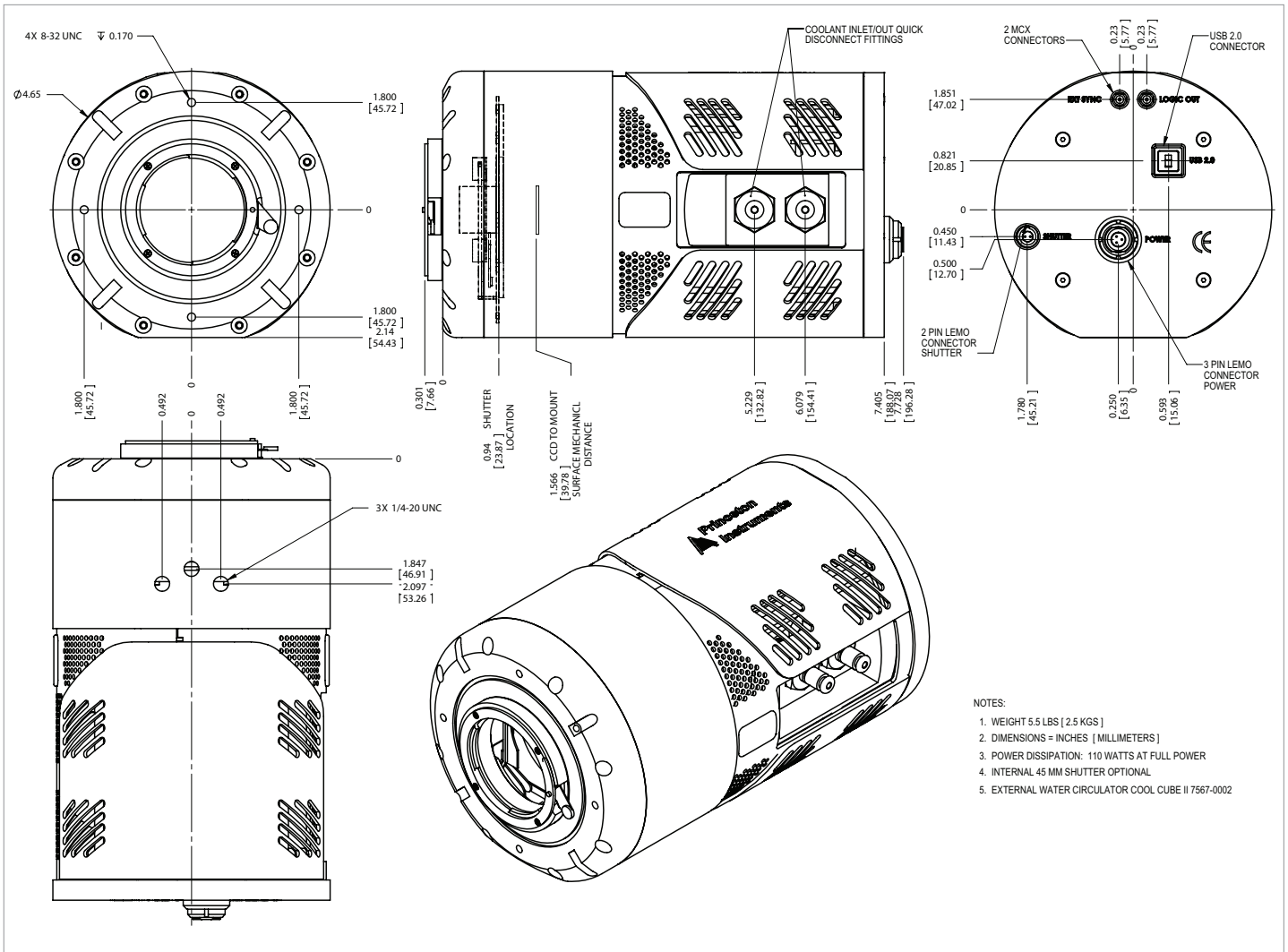

- NOTES:
1. WEIGHT 5.5 LBS [2.5 KGS]
 2. DIMENSIONS = INCHES [MILLIMETERS]
 3. POWER DISSIPATION: 110 WATTS AT FULL POWER
 4. INTERNAL 45 MM SHUTTER OPTIONAL
 5. EXTERNAL WATER CIRCULATOR COOL CUBE II 7567-0002