

COSMOS for Ground-Based Time Domain Astronomy

Introduction

Time domain astronomy is the study of how astronomical objects and unique events **vary over time**. It provides an alternative method to understanding the **extreme phases** of galaxy and stellar evolution, through the investigation of events such as **supernovae**, and gamma-ray bursts, indicative of the collapse of massive stars and the creation of neutron stars or black holes, and objects such as **variable pulsars** and stars, to name a few. Characteristics of these events can be **periodic, random, and bursting in nature**, with variation across all objects and events. This makes time domain astronomy a challenging key field within astronomy and astrophysics.

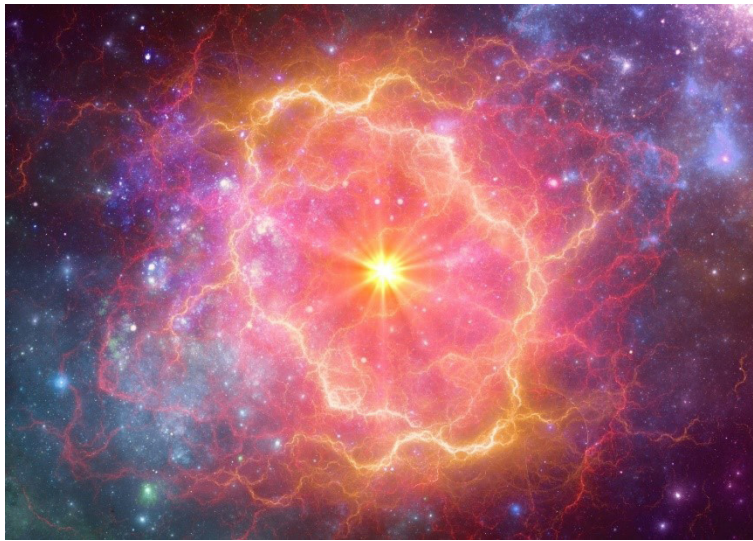


Figure 1: An artist's rendition of a supernovae, an example of an astronomical object of interest for time domain astronomy

Time domain astronomy covers much of the electromagnetic spectrum. It utilizes both **dedicated satellites**, looking at regions of the electromagnetic spectrum that are absorbed by the atmosphere, and **dedicated ground-based**, fully robotic telescopes. These systems work together to capture the optical, infrared, and radio waves, describing the **full electromagnetic profile** associated with each astronomical object or event. Typically, the dedicated satellite will detect a new event, **triggering** the remote, robotic telescope to point at the event (or object) and begin acquiring.

There are different methodologies used for **capturing temporal information** of these variable objects. Imaging is a useful technique for **visualizing** these temporal changes, with spectroscopy typically utilized for looking at the **chemical compositions** of each object or event. Polarimetry is another common technique used within time domain astronomy to understand the **associated variability in magnetic field**. This is particularly useful for gamma-ray bursts, with the polarized light highlighting variations within the magnetic field of the burst.

These events can **vary drastically** in timescale, with some events lasting only a few milliseconds, and others lasting several years. As these events are **unique**, it is essential that the devices used to capture any variation on the very small timescale are optimized.

Camera/Sensor requirements

As there is a lot of variance between these astronomical events, with each being truly unique, the light intensity of the source is **often unknown**. To avoid potentially saturating the sensor, typically a **large number of frames** are taken at short exposure times. Although this method reduces the likelihood of saturation, the shorter exposure time **limits the number of photons** that can be detected. Therefore, a high quantum efficiency is essential to ensure that a high proportion of photons detected are converted into photoelectrons and therefore signal.

A high quantum efficiency is also important as integration over longer periods of time is **not feasible** within time domain astronomy. Once astronomical events are over, they **typically fade**, sometimes within a matter of minutes. A higher quantum efficiency allows for these fainter signals to **still be detected** even though the event has finished (i.e. measurement of the residual signal from the event).

To gain the most out of the information gathered during the fading of an event, the images acquired are typically **stacked**. To maintain a high signal-to-noise ratio of these stacked images, a **low read noise is required**, so that fainter stars or aspects of the event/object can still be detected. When frames are stacked, not only the signal is summed, but **also the read noise**, so low read noise is essential. Computational stacking also allows for atmospheric turbulence to be **partially corrected for**, making a low read noise as integral as a high quantum efficiency.

As these events/objects are unique and imaged on such short timescales, the **true brightness is unknown**. To determine the true brightness, **reference stars**, which have known brightness, are also imaged within the frames. By measuring how the shorter exposure times have influenced the brightness of any reference stars and comparing to the known brightness of the reference star, the true brightness of the event/object can be obtained. The **more reference stars** in each image, the better the **image calibration** can be. Therefore, the larger the physical area of the sensor, the better for image and brightness comparison.

COSMOS for Time Domain Astronomy

COSMOS, the large format, back-illuminated CMOS camera from Teledyne Princeton Instruments, has many of the camera qualities required for time domain astronomy. As it is **back-illuminated**, COSMOS has a high quantum efficiency over the visible range, with **>90% peak quantum efficiency** as shown in Figure 2. This means that it will be able to convert a high percentage of photons into photoelectrons, even with shorter exposure times. COSMOS also has a low read noise, with the ability to achieve **down to 0.7 e- read noise**. This makes the COSMOS suitable for not only imaging the **early stages of an event**, or object, but also detect any **event afterglow**. This low read noise is also advantageous for **stacking of frames**, maintaining a **high signal to noise ratio** even with the summation of frame read noise.

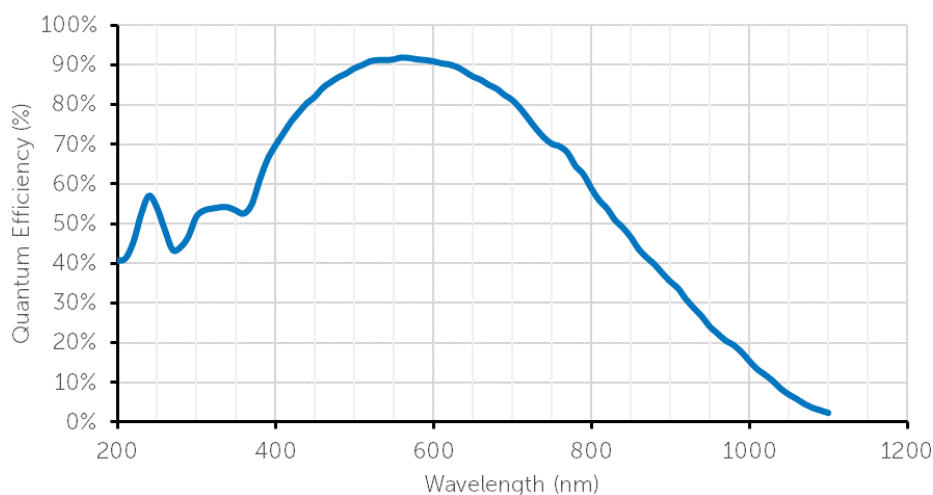


Figure 2: Quantum efficiency curve of the COSMOS camera, showing high quantum efficiency over the visible range and a peak quantum efficiency of >90%.

COSMOS has a **large imaging area**, with 3k x 3k, 6k x 6k, and 8k x 8k formats, all with 10 μm pixels. COSMOS has a sensor diagonal of 43 mm, 92 mm, and 115 mm respectively, as depicted in Figure 3. These large imaging areas allow for the capture of **multiple reference stars** within each frame, ensuring a high level of image comparison and calibration.

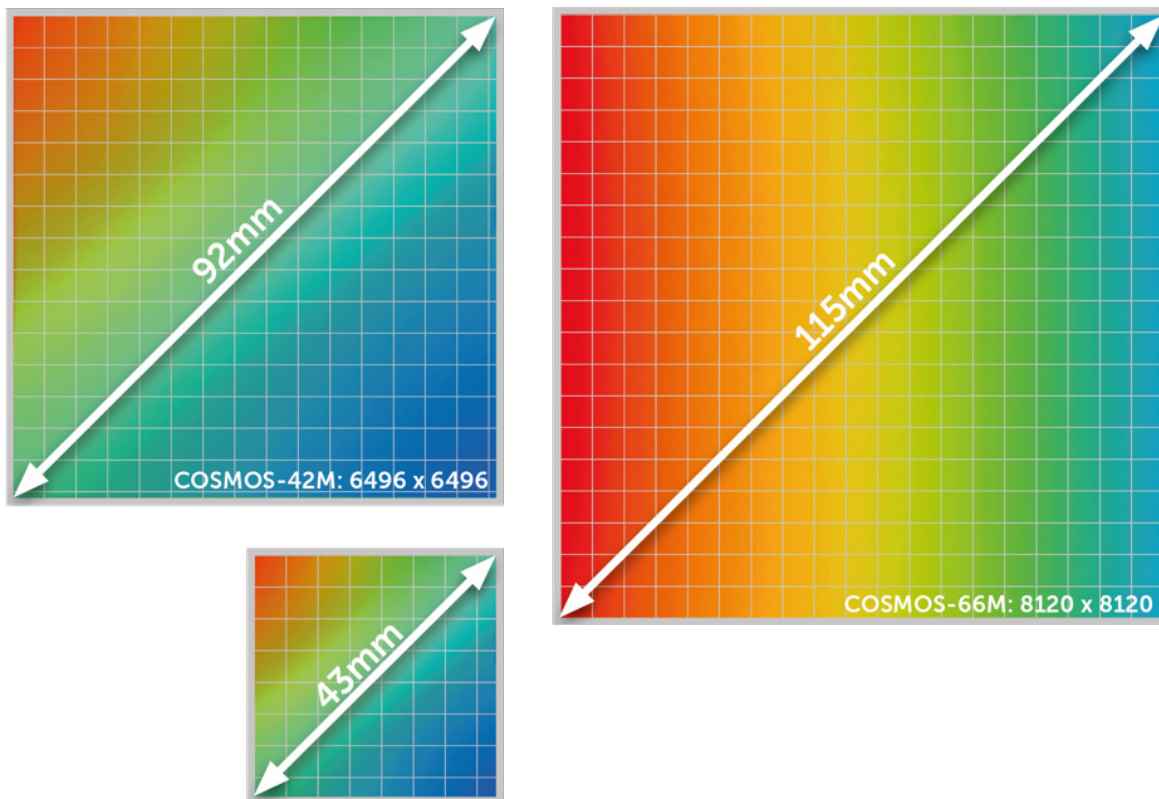


Figure 3: Representative sensor sizes for each of the three COSMOS sensor models, with the 3k x 3k having a sensor diagonal of 43 mm, the 6k x 6k having a sensor diagonal of 92 mm, and the 8k x 8k having a sensor diagonal of 115 mm.

Due to the low dark current and large sensor areas, **full frame CCDs** have been traditionally used within time domain astronomy. In order to completely block any incident light during readout, full frame CCDs utilize a **mechanical shutter**. Mechanical shutters have **finite lifetimes** and often need to be replaced frequently when the camera is in heavy use. This can be problematic for telescopes used within time domain astronomy, as they are typically fully robotic, remote telescopes, meaning that **maintenance can be challenging**.

Additionally, opening and closing a mechanical shutter is relatively slow, introducing **quantitative errors** for the shorter exposure times essential for time domain astronomy. As COSMOS is a CMOS detector, it utilizes a **fast, electronic shutter** where exposure is stopped by shifting detected photoelectrons into a frame storage area before readout. Not only is an electronic shutter **more precise** than a mechanical shutter, but also leads to **less dead** time of the detector, during which the camera is not exposing to light. This means that subsequent exposures can already start as signal is readout from the storage area, allowing **image capture of the entire event**, rather than periodic snippets.

As time domain astronomy relies on the synergy between **dedicated satellites and ground-based telescopes**, it is common for the telescopes to be robotic and remotely controlled. Therefore, it is key that any camera or device can **easily integrate into the existing software**. COSMOS will be completely controlled by PICam, a software development kit from Teledyne Princeton Instruments. **PICam** is available for both **Linux** and 64-bit Windows, allowing complete control of COSMOS in either of these operating systems. PICam offers direct control of the camera, with **flexible configuration and integration** with other languages such as Python. In this way, COSMOS can easily integrate into the software controlling any remote, robotic telescope.

Conclusion

Time domain astronomy is the study of how astronomical objects and events vary and **change with time**. It is a research area that utilizes a lot of **different techniques** to understand the most about each unique event.

Typically, time domain astronomy takes **multiple images with short exposure times** to prevent saturation of the sensor, requiring a camera/device with **high quantum efficiency** and **low read noise** (to maintain high signal to noise ratio when stacking frames). In addition, as each event is **unique**, the true brightness is unknown, so **multiple reference stars** are required to ensure calibration of the object of interest. To measure multiple reference stars within an image, a **large physical sensor area** is required.

The **COSMOS**, with **peak quantum efficiency >90%**, low read noise down to **0.7 e-**, and sensor sizes up to a diagonal of **115 mm at 8k x 8k pixels**, meets the essential parameters for time domain astronomy. In addition, the reduced error from its **electronic shutter**, and **software development kit** for full control with integration into operating systems such as **Linux**, makes the COSMOS optimal for time domain astronomy.